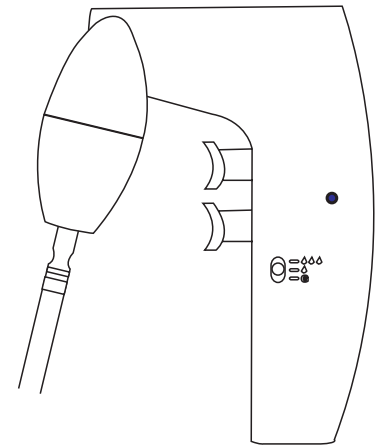
# PROFILLER™ *electro* 446

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#### SAFETY PRECAUTIONS

- Read operating instructions carefully. Observe manufacturer's recommendations.
- Refer to and follow regulations about handling of potentially hazardous reagents.
- Wear protective clothing including goggles and gloves.
- Do not use the Profiller™ *electro* with liquids incompatible with silicone, PP, ABS, ASA and SAN copolymers.
- Do not allow liquid to enter the instrument.
- Never use with highly flammable liquid or in an atmosphere with danger of explosion.
- Should the instrument fail to work to specification, stop pipetting and refer to Trouble Shooting section.
- Before cleaning disconnect from the charger (5).
- Do not clean the main charger.
- All repairs must be carried out by an authorized dealer. Only use the manufacturer's accessories and spare parts.

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#### APPLICATION LIMITS

Profiler™ *electro* 446 is not autoclavable, except the silicone pipette holder (9). Instrument should not be used at temperatures below 10 °C and above 35 °C. This product should be used only for its intended purpose. When instrument is not in use, store on bench stand or wall bracket. Preferably remove wall stand fixture (see Fig. 2) when used on a bench. Charger is only to be used indoors and protected against moisture. Only recharge instrument with manufacturer's original charger.

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#### DESCRIPTION

Profiler™ *electro* 446 is an easy to use and lightweight electronic pipette controller, which facilitates aspiration and distribution of a variety of reagents. The cordless instrument is compatible with glass or plastic pipettes with nominal volumes ranging 1 to 100 mL. The triggers have been designed to eliminate pressure points and control of intake and delivery is therefore made easier. These triggers also control aspiration and delivery speeds.

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## OPERATION

### Recharging Profiller™ *electro* (Fig. 3)

The Profiller™ *electro* operates on two (2 x 1.2V) rechargeable, environmentally friendly NiMH batteries.

- Plug the charger jack (6) into the base of the Profiller™ *electro*.
- Plug the charger (5) into a suitable wall socket. The charger red indicator will light up.
- Allow unit to charge for 14 hours (75 mA/h) from empty to fully charged.
- Once fully charged, the charger automatically switches and the red indicator goes off.

A red battery warning light (4) is triggered whenever the charge is low. Battery should then be recharged.

Pipetting can also take place while the unit is connected to the charger.

### Speed selection:

The Profiller™ *electro* has a conveniently located slide switch (3) which allows the user to select between three modes of operation:

- Switch position **ooo** (top): Aspire and dispense at full speed (blow-out pipette)
- Switch position **o** (middle): Aspire and dispense at reduced speed (blow-out pipette).
- Switch position **G** (bottom): Aspire at reduced speed, dispense by gravity (to deliver pipette).

### Pipette fitting

Carefully attach glass or plastic pipette by inserting it into the holder of the nose cone without using force.

### Aspiration

Hold the instrument in a vertical position with the pipette tip down. Place tip of the pipette into the liquid to be aspirated. Select mode of operation – top, middle for “blow-out” pipette or **G** for “to deliver” pipette. Depress intake trigger (1). The harder the pipetting trigger is pressed the faster the aspiration.

Fill the pipette until the meniscus level has reached the required mark.

### Dispensing

Place pipette against wall of the receiving vessel and:

In “Blow out” mode: Switch (3) on top or middle position, depress the lower trigger (2). The harder the pipetting trigger is pressed the faster the dispense.

In “To deliver” mode: Switch (3) on **G** position, depress the dispense trigger (2). Lower trigger (2) opens the air inlet, allowing liquid to run out by gravity. Wait for the required time indicated on the pipette (i.e. EX + 10 s) for dispensing to be completed.

### Removing the pipette

Hold the pipette and rotate gently while pulling out from the holder.

### TIGHTNESS

Take care of the instrument to keep it in perfect working condition. To test tightness, attach a pipette and fill it at least to half of its nominal capacity. Drops falling from the pipette tip indicate a leak. Unscrew nose cone (10), check pipette holder (9) and membrane filter (8). Reassemble correctly.

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## MAINTENANCE

Before cleaning disconnect from the mains charger (5).  
Do not clean the main charger.

**Do not clean the main charger.**

Do not allow liquid to enter the Profiller™ *electro*. Hydrophobic protection filter (8) would clog and suction be impaired. To replace filter, remove pipette, unscrew nose cone (10) counter clockwise, pull out and replace membrane filter. Reassemble unit. Clean instrument with a damp cloth or with most laboratory disinfectants.

Profiler™ *electro* is not autoclavable, except the silicone pipette holder (9).

## Trouble Shooting

The Profiller™ *electro* is virtually service-free. If the recommendations below do not resolve the problem, please contact your authorized dealer. Do not attempt to open the instrument.

Observation	Possible cause	Action
<b>Aspiration capacity is too slow</b>	Filter blocked, dirty or damaged Batteries low	Clean or replace hydrophobic filter Recharge batteries
<b>Pipette drips</b>	Pipette not properly positioned in pipette holder Pipette holder or pipette damaged	Verify pipette mounting Replace pipette or holder

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## WARRANTY

The Profiller™ *electro* 446 is guaranteed for a one-year period (battery pack 3 months) against any factory or material defects. Damages due to non-respect of manufacturer's instructions are excluded from the warranty. Return the instrument to the dealer from whom it was purchased, after obtaining return authorisation, should regular maintenance not eliminate a detected defect.

**Decontaminate instrument prior to returning it.**

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